


*All enveloped viruses from Annex A, EN14476:2013+A1:2015 including HBV, HCV, HDV, HIV, Coronavirus, Influenza, Rabies, Measles virus.

Oosafe® IVF LABORATORY DISINFECTANTS

Oosafe® DISINFECTANT FOR CO₂ INCUBATORS AND LAMINAR FLOW HOODS

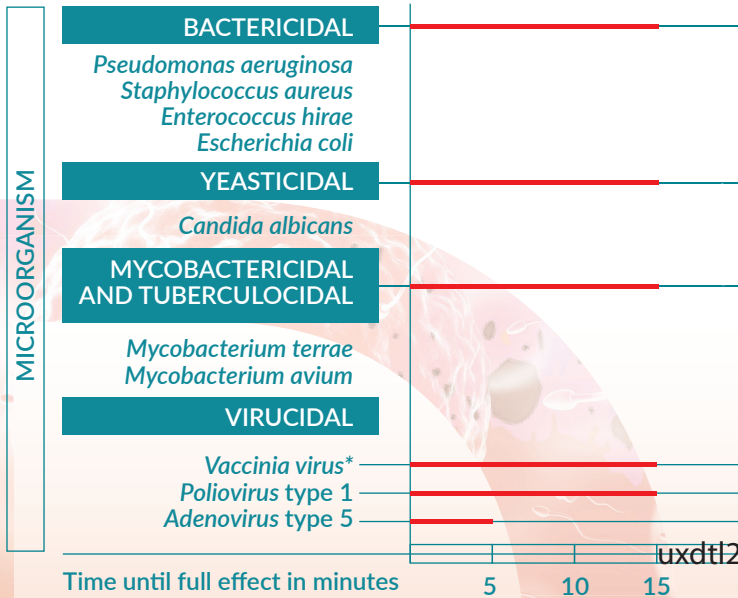

High quality

- Ready-to-use solution
- Anticorrosive
- Securely sealed
- Valid for 3 years, even after opening
- Long term disinfection/low evaporation rate
- Fast disinfection effect
- Transparent and odorless

Health and safety

- MEA and HSSA tested
- Alcohol-free
- No VOC release

Excellent material compatibility

- Stainless steel
- Plastic
- Glass
- Acrylic

Application:

Follow laboratory's surface cleaning procedure before applying the disinfectant.

Spray: Spray and distribute evenly on the surface with a sterile cloth. Wait 15 minutes for the disinfection effect. Wipe the surface to remove any residue.

Wipes: Wipe the surface evenly and wait 15 minutes for the disinfection effect. Close the lid after use. Do not reuse wipes.

Intended use:

Disinfection of CO₂ incubators and laminar flow hoods.



Product reference:
OODIH-0050

Oosafe® Disinfectant for CO₂ Incubators and Laminar Flow Hoods, 50 ml in spray bottle



Product reference:
OODIH-1000

Oosafe® Disinfectant for CO₂ Incubators and Laminar Flow Hoods, 1 Liter bottle with spray



Product reference:
OODIH-5000

Oosafe® Disinfectant for CO₂ Incubators and Laminar Flow Hoods, 5 Liters refill container



Product reference:
OODW-70

Oosafe® Disinfection Wipes, 70 wipes per bucket

Oosafe® IVF LABORATORY DISINFECTANTS

Oosafe® LABORATORY SURFACE DISINFECTANT

High quality

- Securely sealed
- Valid for 3 years, even after opening
- Long term disinfection

Health and safety

- MEA and HSSA tested
- Clear - slightly yellow
- Alcohol-free
- No VOC release

Excellent material compatibility

- Glass, Plastic, Rubber, Stainless steel, Ceramic, Linoleum, Acrylic, Cotton

Features

- Contains non-ionic surfactant to clean and quaternary ammonium compounds to disinfect surfaces at the same time.

Application:

Dilute disinfectant with water according to the table recommendations and apply the diluted cleaning agent on laboratory floors and surfaces. Wait after application for the required disinfection effect and let the surface dry out. If used for laboratory floors, please rinse otherwise it might be slippery.

Intended use:

For disinfection and washing of laboratory surfaces and floors.



Product reference:
ODSF-00050

Oosafe® Laboratory Surface Disinfectant,
50 ml in bottle



Product reference:
ODSF-02000


Oosafe® Laboratory Surface Disinfectant,
2 Liters bottle with measuring cup



Product reference:
ODSF-10000

Oosafe® Laboratory Surface Disinfectant,
10 Liters refill container


*All enveloped viruses from Annex A, EN14476:2013+A1:2015 including HBV, HCV, HDV, HIV, Coronavirus, Influenza, Rabies, Measles virus.

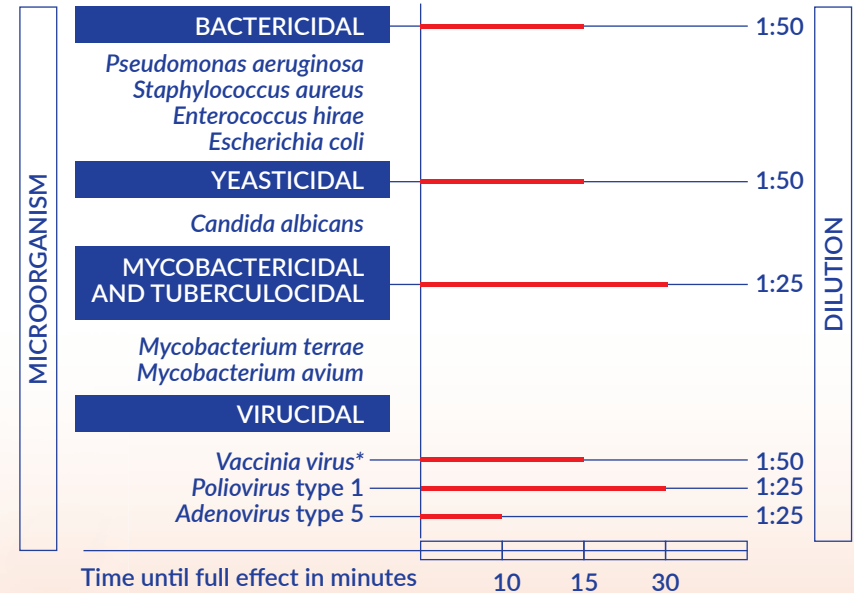

*All enveloped viruses from Annex A, EN14476:2013+A1:2015 including HBV, HCV, HDV, HIV, Coronavirus, Influenza, Rabies, Measles virus.

Oosafe® IVF LABORATORY DISINFECTANTS

Oosafe® HAND DISINFECTANTS

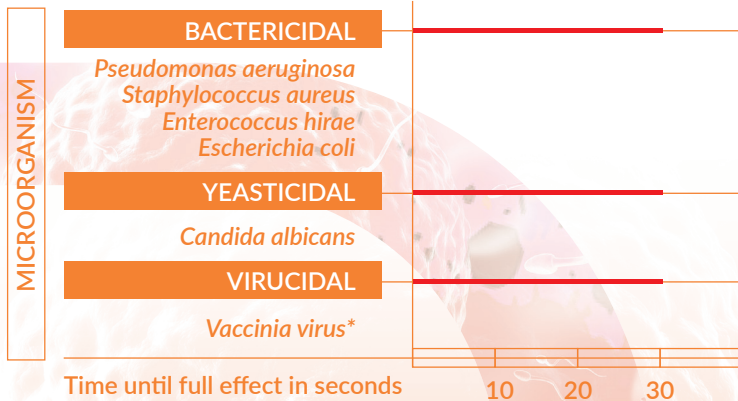

High quality

- Ready-to-use solution
- Securely sealed
- Valid for 3 years, even after opening
- Long term disinfection/low evaporation rate
- Fast disinfection effect
- Transparent and odorless

Health and safety

- Non-sensitizing
- Non-irritating
- MEA and HSSA tested
- Alcohol-free
- No VOC release
- Keeps your skin soft

Application:

Follow laboratory's hygienic procedure for washing hands before using hand disinfectant. Apply minimum 3 ml of disinfectant and rub hands until they are completely dry.

Intended use:

For disinfection of hands.



Product reference:
OODH-0050

Oosafe® Hand Disinfectant,
50 ml in spray bottle



Product reference:
OODH-0500

Oosafe® Hand Disinfectant,
500 ml bottle with pump



Product reference:
OODH-5000

Oosafe® Hand Disinfectant,
5 Liters refill container

ORDER INFORMATION	
Oosafe® Disinfectant for CO₂ Incubators and Laminar Flow Hoods	
Product reference	Product name
OODIH-0050	50 ml in spray bottle
OODIH-1000	1 Liter bottle with spray
OODIH-5000	5 Liters refill container
OODW-70	70 wipes per bucket
Oosafe® Laboratory Surface Disinfectant	
Product reference	Product name
OODSF-00050	50 ml in spray bottle
OODSF-02000	2 Liters bottle with measuring cup
OODSF-10000	10 Liters refill container
Oosafe® Hand Disinfectant	
Product reference	Product name
OODH-0050	50 ml in spray bottle
OODH-0500	500 ml bottle with pump
OODH-5000	5 Liters refill container

